


IR Camera User Guide		
Status: Published		Document ID: 47528
DATE EFFECTIVE 05/18/2021	APPROVED BY Quality Assurance Manager	PAGE 1 OF 2

IR Camera User Guide

1 Setting up the IR Light

- 1.1 Secure the IR light to the moveable arm fixture.
- 1.2 Plug the cable that connects the IR light to the computer into the top of the light. Plug the power cable into the right side of the IR light. See photo below.

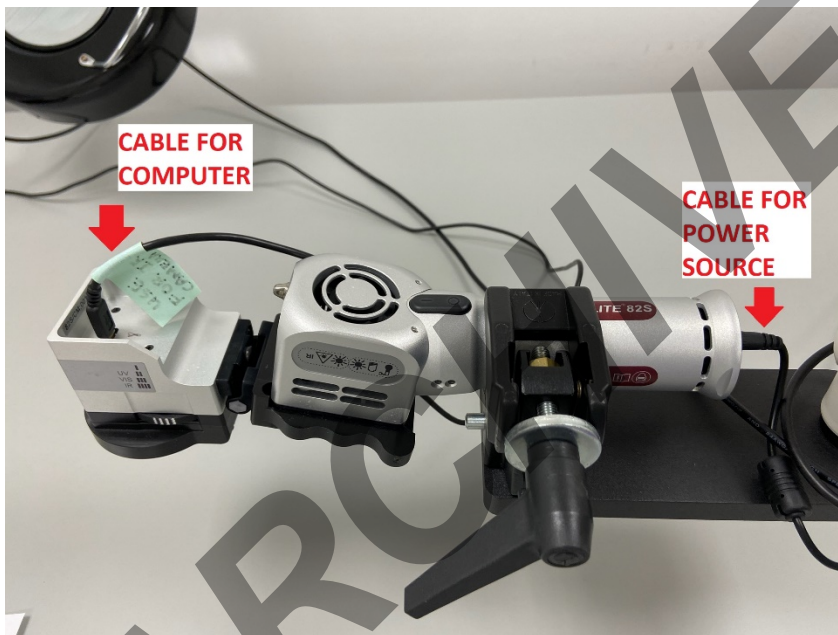

- 1.3 Set the dial on the light to “III”.

2 Using the Crime-Lite Cam Software

- 2.1 In the main window of the software, make sure the “Device Settings” are as follows:

Device: Crime-Lite 82S
Light Source: IR
ML Filter: None
Camera Filter: IR Only III

- 2.2 Set the Exposure setting on the top panel in the software to “Auto”.
- 2.3 To focus the camera:

FORENSIC BIOLOGY EVIDENCE AND CASE MANAGEMENT MANUAL

IR Camera User Guide		
Status: Published		Document ID: 47528
DATE EFFECTIVE 05/18/2021	APPROVED BY Quality Assurance Manager	PAGE 2 OF 2

- 2.3.1 Click “Options” on the top panel in the software.
- 2.3.2 A new panel of settings will appear. Click “Full Screen”.
- 2.3.3 Two settings will appear on the left side of the screen: “Grab Image” and “Auto”.
 - 2.3.3.1 “Grab Image” will take a photo of what is in the camera’s field of view. This photo will appear in the bottom panel of the software, but it does not save anywhere on the computer.
 - 2.3.3.2 Click “Auto” to focus the camera.
- 2.3.4 Deselect “Full Screen” to return to the main window of the software.
- 2.4 Refer to the Evidence Examination manual for how to examine evidence using the IR light.

ARCHIVED