

FORENSIC BIOLOGY PROTOCOLS FOR FORENSIC STR ANALYSIS

Zygem One-Step Sperm Lysis of Sexual Assault Stains or Swabs		
Status: Published		Document ID: 10311
DATE EFFECTIVE 08/04/2017	APPROVED BY Nuclear DNA Technical Leader	PAGE 1 OF 2

Zygem One-Step Sperm Lysis of Sexual Assault Stains or Swabs

Approximately 1/4 of a swab or a 3x3mm cutting of a stain should be used. The Zygem One-Step Sperm Lysis method should only be used for cases involving a male assailant on a female victim. Do not use for male on male cases, female on female cases, nor female assailant on male victim cases.

1 Procedure

- 1.1 Retrieve the following reagents and allow to thaw before use:

Acrosolv
Orange + Buffer
forensicGEM™ enzyme
Ultrapure water

- 1.2 After the reagents have thawed, vortex briefly and spin down for a few seconds.

- 1.3 Create a master mix for n+1 samples (make sure to add two to the number of samples for the two extraction negatives) using the following ratio:

Acrosolv - 10 µL
Orange+ Buffer - 10 µL
forensicGEM™ enzyme - 2 µL
Ultrapure water- 78 µL

- 1.4 The total master mix added per sample is 100 µL. Record the lot numbers.

- 1.5 Retrieve the sample cuttings in 0.2mL tubes. Have a **witness** confirm that the tube labels are correct.

- 1.6 Ensure that all sample cuttings are lightly pushed to the bottom of the tubes (a new pipette tip for each sample can be used for this.)

- 1.7 Add 100 µL of master mix to each sample tube and extraction negative.

- 1.8 Place the tubes into the thermalcycler and close the lid.

- 1.9 Choose the following program on the thermal cycler to extract the samples: Zygem Sperm.

FORENSIC BIOLOGY PROTOCOLS FOR FORENSIC STR ANALYSIS

Zygem One-Step Sperm Lysis of Sexual Assault Stains or Swabs		
Status:Published		Document ID: 10311
DATE EFFECTIVE 08/04/2017	APPROVED BY Nuclear DNA Technical Leader	PAGE 2 OF 2

1.10 The Zygem Sperm program is as follows:

- 52°C for 5 minutes
- 75°C for 3 minutes
- 95°C for 3 minutes
- 4°C for 2 minutes

1.11 Briefly centrifuge the tubes to spin down any condensation.

1.12 The samples can then be aliquoted for Quantifiler Trio.

1.13 Store the extracts at 4°C.

1.14 Reagent Storage:

- 1.14.1 After tubes have been opened, the enzyme and the Acrosolv should be stored at -20°C.
- 1.14.2 The Orange Plus buffer can remain at 4°C for convenience.

ARCHIVED